


ETAERION™

Non-crystalline colloidal nanoparticles


Dynamic light scattering experiments show a single peak with normal distribution*

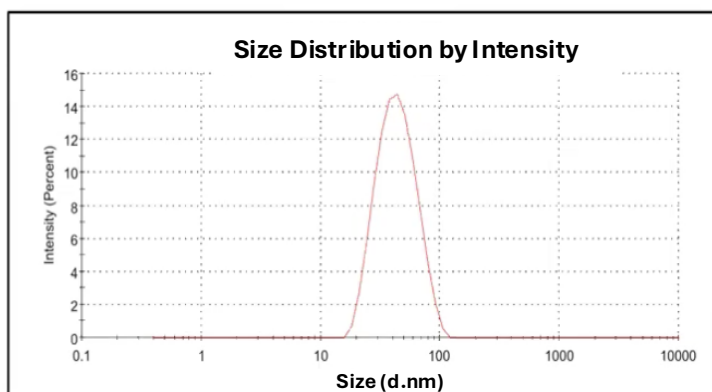


*Crude particles taken directly from the reactor

- Particles are uniform in size and charge
- Particles are not magnetic
- Straightforward to modify post-synthesis

This is a single peak with a negative charge*

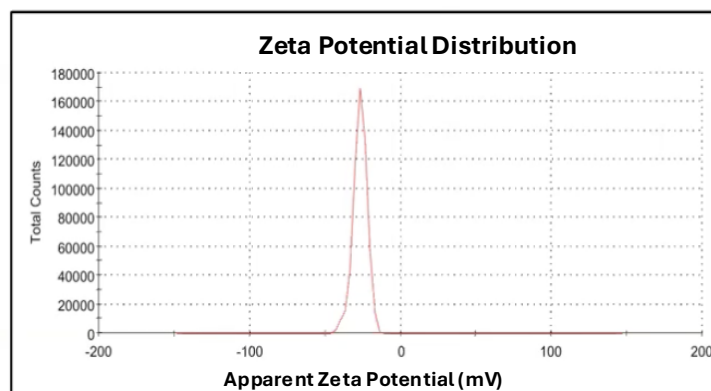

Preservation of ETAERION™ Optical Profile with Dialysis

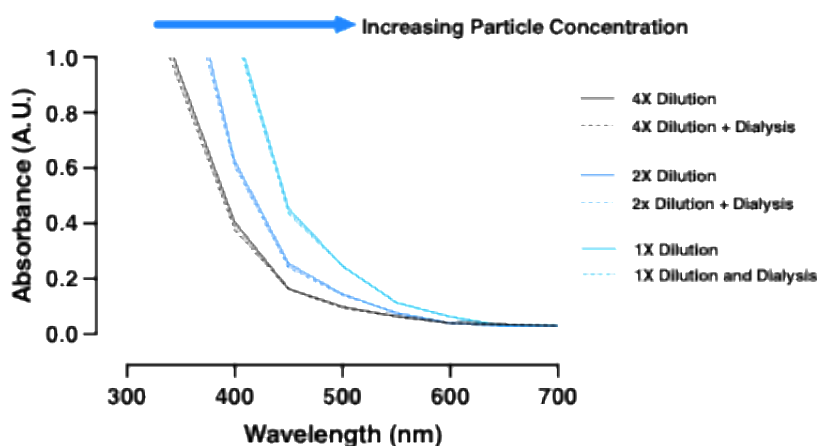

Interpretation: ETAERION™ remains structurally intact and does not release diffusible iron under buffer exchange conditions.


Dialysis against PBS does not alter the absorbance profile, indicating the signal is dominated by particle-associated iron and diffusible iron species are minimal under these conditions. Data suggest ETAERION™ remains structurally intact and does not release diffusible iron under buffer exchange conditions. Data also suggest there are minimal free iron residuals present post synthesis. Absorbance scales with particle concentration.

ETAERION™

Mineral Nanoparticle Colloid

ETAERION™ improves how cells respond to high acute stress


ETAERION Particles Enhance Resilience in Stressed HEK293T Cells

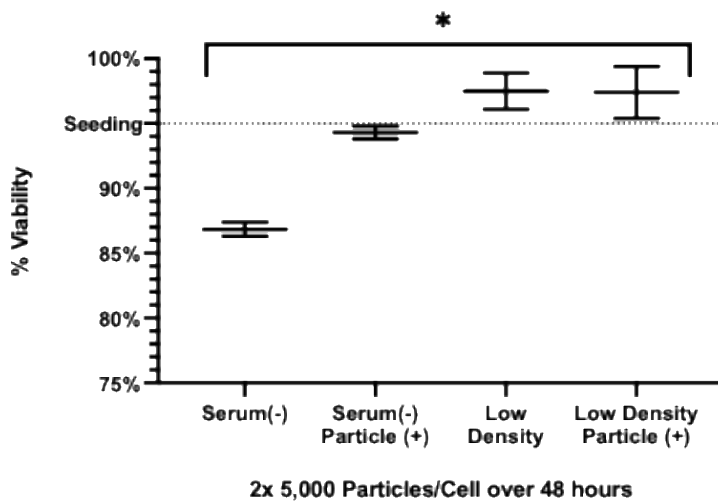

- When serum-starved HEK293 cells are treated with 5,000 particles/cell (every 24 hours for 48 hours,) viability increases
- ETAERION™ particles do not appear to have an effect on cells seeded at low density, suggesting that increased viability is affected by the nature/severity of cell stress

An ordinary one-way ANOVA was performed to compare HEK293T cell viability across four treatment conditions (n = 8 total). A significant difference among group means was observed ($F(3,4) = 15.36, p = 0.0116$), indicating that treatment condition significantly affected cell viability. The model explained a large proportion of the variance ($R^2 = 0.92$). Variance across groups was not homogeneous, as indicated by a significant Brown-Forsythe test ($p < 0.0001$), suggesting that treatment effects differ not only in mean viability but also in variability. These results support a strong, condition-dependent effect of ETAERION exposure on cell viability under stress, consistent with a non-uniform, state-dependent biological response.

Interpretation: ETAERION™ improves viability selectively under stress conditions rather than acting as a universal enhance

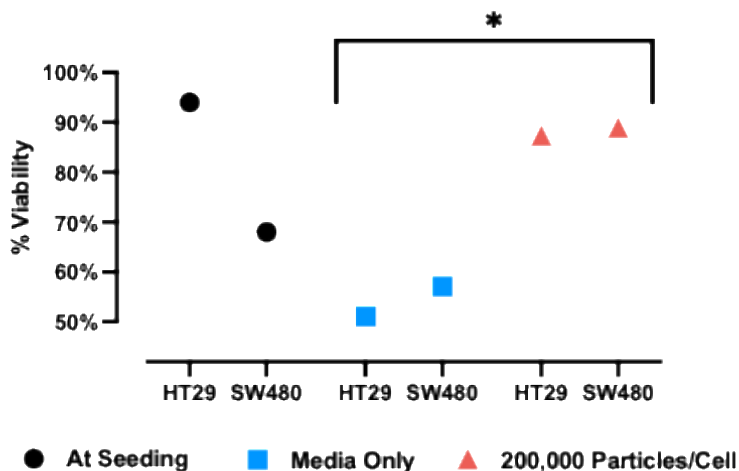


- When HT29 and SW480 cells are treated with 200,000 particles/cell (every 24 hours for 48 hours) viability increases
- Data confirm preservation of viability without driving cell growth across multiple cell lines
- Viability is shown at seeding (black circles), after 48 hours in with media alone (blue squares), after 48 hours with ETAERION™ (red triangles)

Paired two-tailed t-test comparing C002-treated cells to media-only controls across matched colorectal cancer cell lines (HT29, SW480). C002 exposure resulted in a significant increase in viability (mean difference = +34%, $p = 0.041$ (exploratory analysis, n = 2 paired lines).

ETAERION™

Mineral Nanoparticle Colloid


ETAERION™ reduces media acidification in HEK293 cells in acute settings (48 hours)

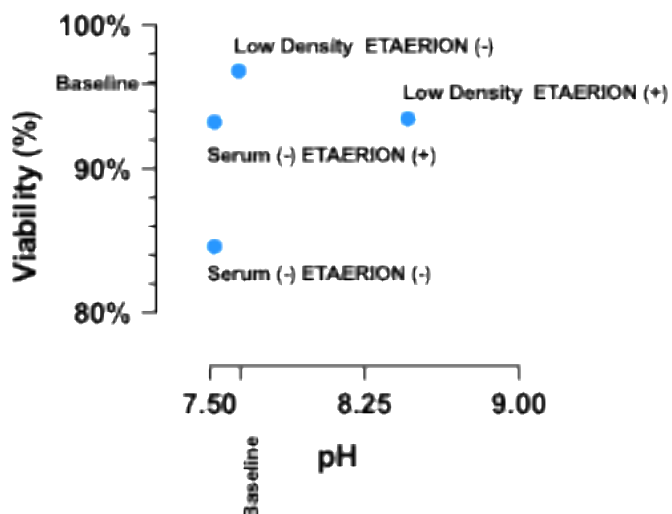

- Serum-starved cells show increased viability when exposed to ETAERION™ (5,000 particles per cell over 48 hours)
 - The magnitude of the effect is reduced relative to prior experiments, consistent with lower particle dose and higher baseline viability
- In low-density conditions, ETAERION™ does not significantly alter viability but increases media pH, indicating reduced acidification

Interpretation: ETAERION™ modulates extracellular pH, with downstream effects on viability dependent on stress context (impact is greater for low-stress conditions; i.e., low-density plating)

Recoverability of HEK293T Cells Under Differing Chronic Stress Conditions

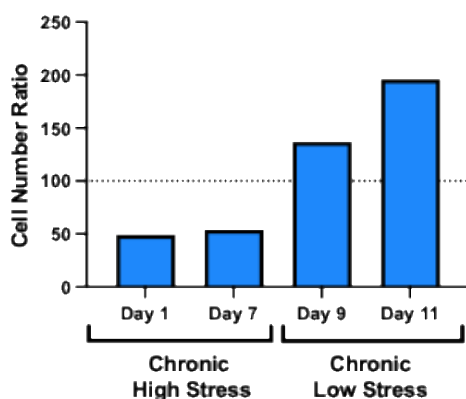

- HEK293T cells with or without 5,000 particles/cell of ETAERION™ were serum starved for seven days. Viability at each time point is shown as a ratio of the number viable cells in the ETAERION™ cultures to the number of viable cells in the serum-only cultures
 - Viability on Day 1 and Day 7 suggest that ETAERION™ confers a disadvantage to cells under chronic high-stress conditions
- On day 7, cells were passed and supplied with normal media
 - At Days 9 and 11, restoration of viability is enhanced for ETAERION™-containing cultures. This crossover behavior is consistent with a state-dependent response


Interpretation: ETAERION™ does not universally enhance growth, but instead modulates system performance within a defined window of viability, where it can enhance recovery following a transition out of chronic high stress environments

ETAERION™

Mineral Nanoparticle Colloid

Under Chronic Stress Conditions (96 hours) there is a bifurcated, cell-stress threshold-based response to EATERION™

In Chronic Setting There is a Bifurcation in Resonse to 5K ETAEION Particles per Cell

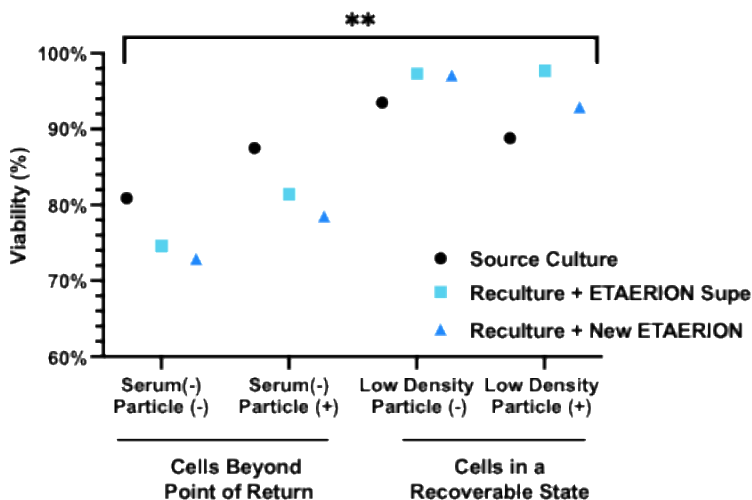

Two-way ANOVA (ordinary, $\alpha = 0.05$) was used to assess the effects of culture condition (row factor) and treatment (column factor) on cell viability. A significant main effect of culture condition was observed ($F(3,6) = 13.79$, $p = 0.0042$), accounting for ~86% of total variance, indicating that baseline stress state strongly determines viability. No significant main effect of treatment alone was detected ($F(2,6) = 0.40$, $p = 0.687$), suggesting ETAERION does not uniformly alter viability across all conditions. Residual variance was low ($MS = 18.6$), consistent with a structured, condition-dependent response. These data support a state-dependent biological effect, in which ETAERION's impact emerges only within specific stress regimes rather than as a global viability enhancer.

- A source culture of HK293 cells was grown for 48 hours in one of four conditions (either serum-starved or low-density conditions with or without particle). Black circles show viability at 48 hours
 - At 48 hours, cells in each condition were re-cultured with either new ETAERION™ (square) or supernatant from the low-density + particle source culture (triangles)
 - When cells were re-cultured, stress conditions (serum(-) or low-density plating) were maintained
- 48 hours following initiation of re-culture, viability was measured
 - In serum-starved conditions, ETAERION™ reduces viability, indicating a reversal of effect under more severe stress
 - ETAERION™ increases viability in low-density (recoverable) cultures
- Response to conditioned supernatant suggests involvement of soluble or environment-mediated effects

Interpretation: These results support a model in which ETAERION™ modulates cellular outcomes in a context-dependent manner rather than as a universal viability enhancer. In addition, results suggest a physicochemical environmental effect may contribute to ETAERION™ activity.

ETAERION™

Mineral Nanoparticle Colloid

ETAERION™ Improves recoverability of cells following freeze-thaw

- Pre-conditioning cells with ETAERION™ 48 hours prior to freezing improves post-thaw recoverability relative to untreated controls
- Inclusion of ETAERION™ in freeze media does not improve outcomes and may transiently reduce viability immediately after thaw
- Addition of ETAERION™ post-thaw enhances recovery, with the greatest benefit observed in cells pre-conditioned prior to freezing
- The strongest recoverability is observed when ETAERION™ is applied both before freezing and after thaw, indicating a staged effect on cellular resilience and recovery